

Preparing M2

1.0 Equipment

1.1 Fridge

2.0 Supplies

2.1 M2

2.2 14ml falcon tubes

2.3 0.2 μ m syringe filters

2.4 Luer Lock 20ml syringe

3.0 Procedure

3.1 Bottles of 50ml M2 medium are stored in fridge until required. Check expiry date displayed on bottle before use.

3.2 Label 4 x 14ml falcon tubes with 'M2' and the date.

3.3 Filter M2 using a 20ml syringe and 0.2 μ M filter.

3.4 M2 can be stored in fridge 5 for up to 7 days.

